

ProbeSure™ COVID-19 One Step RT-PCR Singleplex Kit

Introduction

The ProbeSure™ COVID-19 One Step RT-PCR Singleplex Kit is for the qualitative detection of RNA of the novel Coronavirus (SARS-CoV-2) in human respiratory samples. The kit contains assays for two target sequences in the nucleocapsid (N) gene region and an internal RNase P control assay. The results can be analysed in endpoint or real-time. For full details, please refer to the ProbeSure™ COVID-19 One Step RT-PCR Singleplex Kit User Guide at www.3crbio.com.

Included in the kit

- ProbeSure™ one step RT-PCR Master Mix (2x)
- 20x Assay Mix (Blue lid) for N Gene targets N1 (reports with FAM)
- 20x Assay Mix (Yellow lid) for N Gene targets N2 (reports with FAM)
- 20x Assay Mix (Orange lid) for RNase P genes (reports with FAM)

Not included in the kit

- PCR plate and optically clear seal
- PCR-grade water
- Template RNA and +/- Controls

Storage

All components must be stored at -20 °C upon arrival. Avoid repeated freeze / thaw cycles.


Safety warnings and precautions

This product should only be handled by trained laboratory personnel. Wear suitable PPE when using the product. In case of contact with skin or eyes, wash immediately with water.

	96-well plate	384-well plate
2x ProbeSure-RT Master Mix	10.0	2.5
20X Assay Mix for N1 or N2 or RNase P ¹	1	0.25
Water	<i>Variable</i>	<i>Variable</i>
RNA ²	<i>Up to 9.0</i>	<i>Up to 2.25</i>
TOTAL	20.0	5.0

Table 1. Reagent volume for assembly of reaction mix. Total volumes indicated are recommended for the associated PCR plate type. Include positive, negative, and no-template controls. ¹ A separate total reaction mix must be assembled for each of the three assays: N1, N2 and RNase P. ² Add the maximum volume of RNA possible to maximise RNA copy number in the reaction. However, this should be determined experimentally as some crude extraction methods may introduce PCR inhibitors.

ProbeSure™ Covid-19 process flow

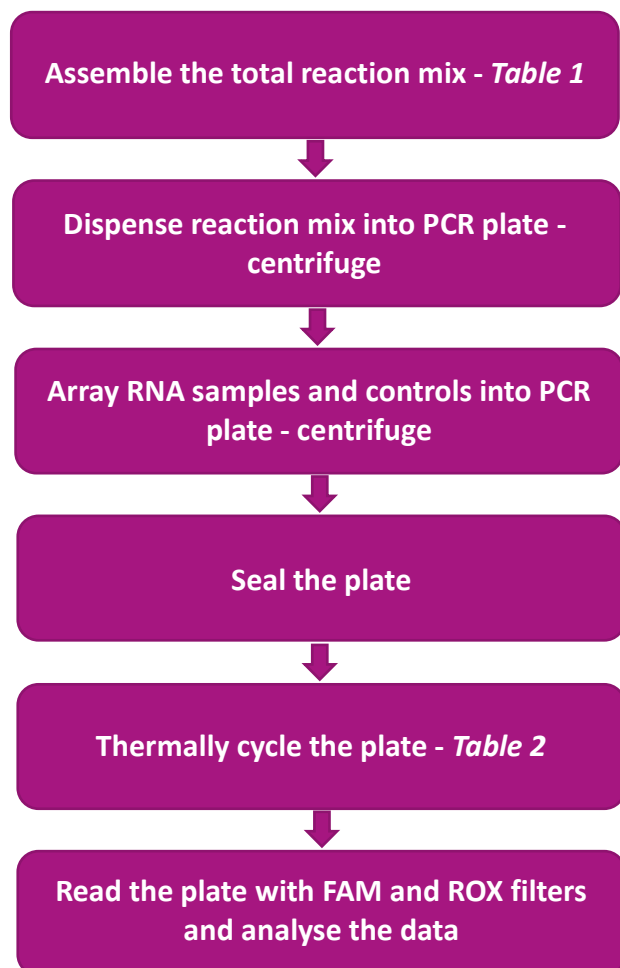

Step	Description	Temperature	Time	N°. Cycles
1	Reverse Transcription	50°C	10 min	1
2	Enzyme activation	95°C	10 min	1
3	Template denaturation	95°C	15 secs	50
	Annealing and extension	60°C	60 secs	

Table 2. Thermal cycling conditions for ProbeSure™ COVID-19 one step RT-PCR Singleplex Kit. PCR can be carried out on any Peltier or water bath thermal cyclers.

ProbeSure™ COVID-19 One Step RT-PCR Singleplex Kit

Data analysis

After thermal cycling, read the FAM and ROX fluorescence in a FRET-capable plate reader. Data can be analysed in real-time or endpoint.

If analysing in endpoint, import the data into a genotype cluster analysis software package. Using the Singleplex kit, the data must be viewed as three separate cluster plots: one for N1 (FAM), one for N2 (FAM) and one for RNase P (FAM) control. For a sample to be COVID-19 positive it would need to be positive on the N1, N2 and RNase P data plots, see *Table 3*.

N1	N2	RNase P	Result
+	+	+	COVID-19 positive
-	+	+	COVID-19 negative
+	-	+	COVID-19 negative
-	-	+	COVID-19 negative
-	-	-	Reaction failed
+	+	-	False positive
+	-	-	False positive
-	+	-	False positive

Table 3. Analysis of the results of the cluster plots of the three singleplex assays of the ProbeSure™ COVID-19 one step RT-PCR kit.

Ordering information

Please visit www.3crbio.com

Licence information

For Research Use Only. Not for use in diagnostic procedures.

3CR Bioscience Ltd disclaims all warranties with respect to this documentation. The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly or by implication. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration.

©2020 3CR Bioscience Ltd. All rights reserved. Intended for molecular biology applications.

For any queries about this guide please contact:

Email: support@3crbio.com

Tel: 01279 940 983

www.3crbioscience.com